NETTOYEUR A ULTRASONS
REF 1058453

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ULTRASONS CLEANER

OPERATION

Elmasonic S350 R "Lab Technology" allows the cleaning of test sieves, degassing HPLC solvents and the processing of analytical samples using state-of-the-art ultrasound technology.

Equipped with microprocessor-driven ultrasonic modes and various specifically defined programs, this device is a real support in carrying out the daily tasks of laboratory cleaning.

The proper functioning of a sieve analysis depends beforehand on the impeccable cleanliness of the mesh and involves the total elimination of all grains or residues.

How to clean the sieve: insert the test sieve with its support into the tank.

The special "sieve cleaning" function allows, thanks to the constant switching between the different ultrasonic modes, to generate strong pulses and thus obtain optimum cleaning results.

HPLC solvents are degassed in Erlenmeyer vials or solvent-resistant 1-litre bottles.

At the back of the Elmasonic S350R is a fastener to attach a laboratory tripod.

The container thus attached will be immersed in the tank previously filled with water and detergent.

The special "Degas" function will allow for about 30 minutes, to degass the solution effectively.

During this process, the gas first forms large macroscopic bubbles, which will be expelled on the surface during short pauses in sound activity.

The "Degas" mode is also particularly well suited to sample degassing, e.g. for carbonic acid degassing in the food sector.

The additional "sample prep" function is intended to perform ordinary tasks such as mixing, dissolving, scattering, cleaning etc.

Technical characteristics

Automatic programs for sieve cleaning and processing test samples

Effective and silent "Degas" function (Gas-Liftout effect)

37 kHz ultrasonic transducers

Cavitation-resistant stainless steel tank with filling level marker (3, 4, 5 L)

Useful space for sieve up to 500 mm

Short-term operation for 1 to 30 minutes, or continuously, adjustable by rotary switch

LED ad for deposit value and remaining cleaning time

Connection for fixing to a tripod (silent operation)

Detachable electric cord equipped with a safety plug

Technical data

Mains voltage (Vac) 1	15-120 V / 220-240 V	Max. filling capacity (liter.) 57.0
Ultrasonic frequency (kHz)	37	Weight (kg) 21.0
Total absorbed power (W)	600	Stainless steel material
Effective ultrasonic power (W)	600	Stainless steel housing material
Maximum peak ultrasonic power * (W)	2400	Handles ✓
External dimensions of device W / H (mm)	650 / 455	CE compliant ✓
Internal tank dimensions W / H (mm)	600/245	Protection class IP 20
Internal basket dimensions W / H (mm)	545/70	
Useful space Ø for sieve up to (mm)	500	

** The choice of waveform has been adapted to the size of the tanks.*

Due to the shape of the waves, factor 4 results for the maximum peak value of the ultrasonic power.

Additional accessories - contact us:

Bucket **ref 106 2233**

Plastic cover **ref 106 3389**

Support for test sieve - **on request**

Additional accessories **on request**